

# NEWSLETTER

**ALFA WASSERMANN**  
Separation Technologies, LLC

Issue 4 | August | Year 2025

## New tool announcement – Holder, Top Rotor Nut (P/N 506360)

To simplify your installation process, we're introducing a new helper: the Holder, Top Rotor Nut, now included with all new units.

This lightweight, easy-to-use tool keeps the top nut in place during shaft installation — saving time, freeing your hands, and improving precision.

### The problem:

Traditionally, installing the top shaft onto the top cap required holding the nut in place by hand — not only awkward, but it also limited visibility and control.



The installer must hold the nut manually while guiding the shaft into position.

### The solution:

Our new Holder, Top Rotor Nut (P/N 506360) solves this. It clips securely around the nut, holding it in place during installation — no hands needed. With the holder installed, the nut stays perfectly aligned — giving you full visibility and both hands free.



A compact, purpose-built part to hold the rotor nut in position.

### Key Benefits

- No need to use your fingers to hold the nut
- Frees both hands for accurate shaft alignment
- Reduces risk of misalignment or cross-threading

### Interested in the Holder, Top Rotor Nut?

Please contact our sales department at:

- [bvestal@alfawassermannus.com](mailto:bvestal@alfawassermannus.com)
- [rfreijer@alfawassermann.nl](mailto:rfreijer@alfawassermann.nl)



[www.awst.com](http://www.awst.com)

This document contains technical information and data proprietary to Alfa Wassermann Separations Technologies. The disclosure to others or use of this technical information or data in any way without prior written consent from Alfa Wassermann Separation Technologies, is hereby expressly prohibited.

Page 1 of 1

Units involved: ☒ Air-(p)KII ☒ e-(p)KII ☒ Promatix ☐ AFH